

Procédure de prélèvement




Placer le gel au congélateur au moins 48 h avant l'expédition



1. Mettre les gants




2. Délimiter la zone à tester

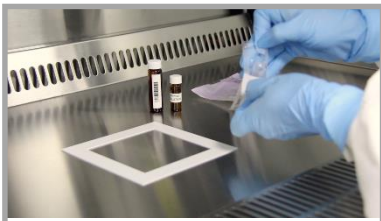


3. Déballer la pipette stérile




4. Déballer les écouvillons stériles

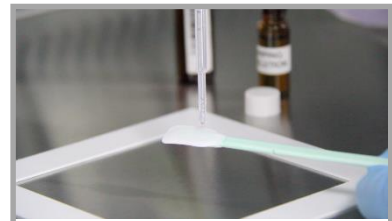

5. Humecter l'écouvillon avec 3 gouttes de solution

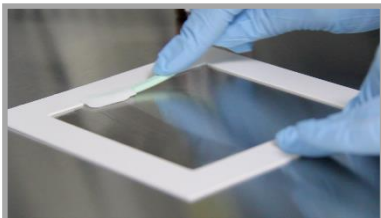

6. Essuyer la surface verticalement

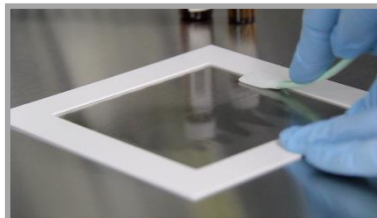

7. Tourner l'écouvillon et essuyer la surface horizontalement




8. Insérer l'écouvillon dans un tube préalablement ouvert et casser le manche



9. Refermer le tube



Prélèvements de surfaces non planes possibles

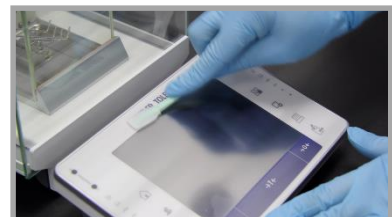

Prélèvements sur des écrans tactiles possibles

Répéter les étapes 2 à 9 pour les autres prélèvements

Consulter la vidéo sur cytoxlab.ch/infos-pratiques

Wipe sampling procedure




Put the ice pack in the freezer
48 h before shipping



1. Put on the gloves




2. Define the surface to be
tested

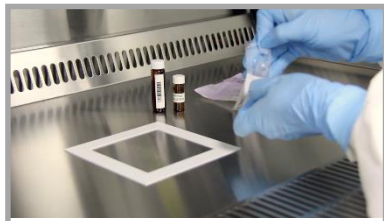


3. Unpack the sterile pipette




4. Unpack the sterile swabs

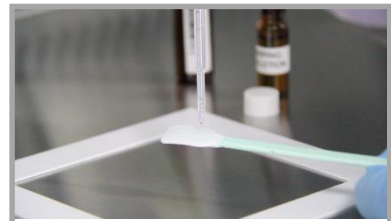

5. Moisten a swab with 3
drops of sampling solution

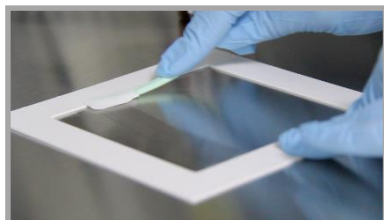

6. Wipe the surface vertically

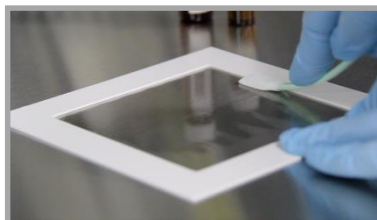

7. Turn the swab and wipe the
surface horizontally




8. Insert the swab in the
previously opened tube and
break the handle



9. Close the sample tube



Wipe sampling of non-planar
surfaces possible

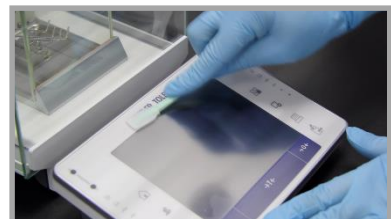

Wipe sampling of touch
screens possible

Repeat the steps from 2 to 9 for additional samples

Consult the video on cytoxlab.ch/infos-pratiques